


## Smart Cart Charging Garage

ME-1243

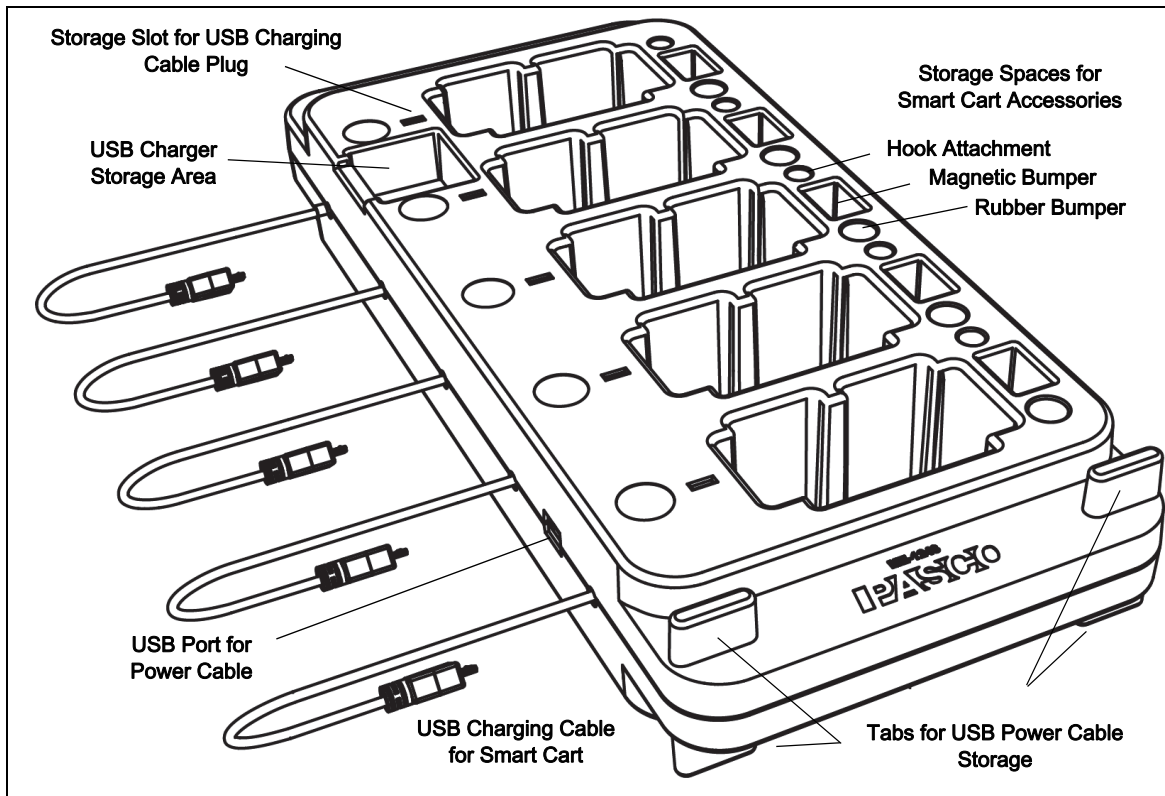

### Introduction

The PASCO ME-1243 Smart Cart Charging Garage is designed to work with the PASCO ME-1240 or ME-1241 Smart Carts. The Charging Garage can hold up to five Smart Cart's at one time. It takes approximately five hours to fully charge five Smart Carts.

#### PS-2575 Single Port USB Wall Charger

The Smart Cart Charging Garage includes a two meter long USB cable (USB-A to Micro-B) and the PS-2575 Sin-

gle Port USB Wall Charger. The charger includes a power adapter and four interchangeable plugs.



The Smart Cart Charging Garage has built in storage spaces for the Smart Cart Accessories (Hook Attachment, Magnetic Bumper, and Rubber Bumper). The garage also

has a storage area for the single port USB wall charger. Along the top of the garage are five small slots for storing the charging plugs when they are not being used.



The top illustration shows four Smart Carts in position with their USB Charge Cables attached. The fifth charger cable is stored in the slot on top of the garage. The USB Power Cord is shown connected to the USB Port on the side of the Charging Garage.

The bottom illustration shows the storage places for the Smart Cart Accessories.

## Operation

Connect the larger end of the USB Power Cord to the USB port on the end of the power adapter. Connect the smaller end of the USB Power Cord to the USB Port on the side of the Smart Cart Charging Garage. Plug the power adapter into a compatible electrical receptacle.

On the Smart Cart, push the plunger all the way into the body of the cart before putting the cart into one of the slots of the Charging Garage. Connect the USB Charging Cable into the USB port on the side of the Smart Cart. The battery status LED on the Smart Cart will indicate what is happening to the battery. The battery status shines yellow when the battery is charging, and shines green when the cart is

charged. Insert as many as five Smart Carts. The Smart Cart Charging Garage has a fixed regulated power supply so it does not matter how many of the five slots are filled with Smart Carts.

The Smart Cart's charging circuitry will disconnect itself from the battery when fully charged.

## Storage

Smart Carts can be stored in the Smart Cart Charging Garage for long periods of time without harming the battery. The Smart Cart circuitry monitors the battery level, and when it falls below a certain level, the circuitry will start charging the unit. When the battery is fully charged, the Smart Cart charging circuitry will stop charging the unit.



## Technical Support

For assistance with any PASCO product, contact PASCO at:

Address: PASCO scientific  
10101 Foothills Blvd.  
Roseville, CA 95747-7100

Phone: 916-462-8384 (worldwide)  
800-772-8700 (U.S.)

Web: [www.pasco.com](http://www.pasco.com)

Email: [support@pasco.com](mailto:support@pasco.com)

## Product End of Life Disposal Instructions:

This electronic product is subject to disposal and recycling regulations that vary by country and region. It is your responsibility to recycle your electronic equipment per your local environmental laws and regulations to ensure that it will be recycled in a manner that protects human health and the environment. To find out where you can drop off your waste equipment for recycling, please contact your local waste recycle/disposal service, or the place where you purchased the product.

The European Union WEEE (Waste Electronic and Electrical Equipment) symbol (to the right) and on the product or its packaging indicates that this product **must not** be disposed of in a standard waste container.

## More Information

For more information about the latest revision of this Instruction Sheet, visit:

[www.pasco.com/manuals](http://www.pasco.com/manuals)

and enter the Product Number.

For information about the Smart Cart Charging Garage or any PASCO product, what software to use, and what PASCO tracks and other accessories are available, check the PASCO web site:

[www.pasco.com](http://www.pasco.com)

## Warranty, Copyright, and Trademarks

**Limited Warranty** For a description of the product warranty, see the PASCO catalog. **Copyright** The PASCO scientific 012-15277A *Smart Cart Charging Garage Instruction Sheet* is copyrighted with all rights reserved. Permission is granted to non-profit educational institutions for reproduction of any part of this manual, providing the reproductions are used only in their laboratories and classrooms, and are not sold for profit. Reproduction under any other circumstances, without the written consent of PASCO scientific, is prohibited. **Trademarks** PASCO and PASCO scientific are trademarks or registered trademarks of PASCO scientific, in the United States and/or in other countries. All other brands, products, or service names are or may be trademarks or service marks of, and are used to identify, products or services of, their respective owners. For more information visit [www.pasco.com/legal](http://www.pasco.com/legal).